

BT100S Variable Speed Stainless Steel Peristaltic Pump Serials



Product Description

BT100S Variable Speed Stainless Steel Peristaltic Pump Serials

BT100S speed variable peristaltic pumps adopts imported core processor, high quality stepper servo motor drive, speed control accuracy is $\pm 0.2\%$. Through the APP software, the pump can be remotely controlled and the running status can be monitored in realtime. RS485 communication, Modbus protocol, can be connected to the upper computer such as computer, single chip microcomputer, PLC, frequency converter, man-machine interface through a variety of signal modes, to achieve systematic centralized control.

Applications:

Liquid addition for synthetic reactions, Liquid discharge, TFF ultrafiltration, Liposuction, Coating and granulation (for seeds and chemical/pharmaceutical applications), Supporting fermentation equipment, Chromatography columns, Laboratory

reaction vessels, Liquid circulation, Supporting filling machines, Coating (for fiberglass electrostatic agents, enamel wire lubricants, and handheld device adhesives)

Features:

1. LF-LCD-OS software system, high definition lattice LCD display, switching between Chinese-English interface, simple parameter setting, intuitive state display.
2. High-quality and high-torque stepper servo motor drive, speed accuracy, wide range, operation stability, high-precision flow transmission.
3. Industrial mask key operation, simple and convenient, high durability.
4. LF-Easy Change universal design, large torque output, strong expansibility, can match various high-performance pump head.
5. Start/Stop, adjust speed, reversible direction, full speed and, state memory (power-down memory).
6. The parameters of running time, interval time and cycle times can be set to meet the requirements of timing, quantitative, liquid dispensing and flow testing.
7. Slow speed stop and suction function, which can effectively prevent the dropping of liquid when the machine stops.
8. Keyboard lock function to prevent misoperation.
9. The remote start-stop, adjust speed and timing operation can be realized by using Lead Fluid APP software. It also has monitoring functions such as stop alarm, change of pump tube and so on (optional).
10. Stainless steel 304 material housing, effective prevention of erosion of various organic solvents, easy to clean.
11. The internal structure adopts double-layer isolation design, and the circuit system provides special protection, which effectively improves the anti-dust, moisture-proof, anti-corrosion and oxidation resistance.
12. Super anti-interference characteristic, wide voltage design, suitable for complex power supply environment.
13. RS485 communication interface, Modbus protocol, easy to connect with various control devices.
14. Acceptable multiple external industrial control signals to control start-stop, reversible direction and speed adjustment, optically coupled isolator.
15. Support start-stop, reversible direction, current speed etc. working state signal output.
16. External thermal printer (optional), real-time printing operation parameters.

17. It can be equipped with foot switch, pulse damper, leak detector and other accessories.

Specifications:

1. Flow range: 0.0001~750 mL/min
2. Speed range: 0.1~150rpm
3. Speed resolution: 0.1rpm (speed≤100rpm), 1rpm (speed>100rpm)
4. Speed accuracy: <±0.2%
5. Timing function: Timing running time 0.1~999 S/Min/H/D, interval time 0.1~999 S/Min/H/D
6. Interval time: 0.1~999.9 S/Min/H, time unit adjustable
7. Cycle times: 1~999, “0” Infinite cycle
8. External control signal input: (1) Start/Stop: passive contact, external control input level (5~24V); (2) Forward/Reverse: passive contact, external control input level (5~24V); (3) Speed Regulation: analog 0~5V, 0~10V, 4~20mA can be set
9. External control signal output: (1) Start/Stop: Level Signal (following input voltage) (2) Forward/Reverse: Level Signal (following input voltage)
10. Communication interface: RS485 communication interface, Modbus protocol is available, WiFi function (optional)
11. Power supply: AC 100~240V, 50Hz/60Hz
12. Power consumption: <40W

Order Number	Channels	Pump Head	Tube Size	Flow Rate of Single Channel (ml/min)
BT100S-D1	1	DG6-1/DS6-1 (6 Rollers)	W.T.:0.8~1mm, I.D.:≤3.17mm	0.00016~49
BT100S-D2	1	DG10-1/DS10-1 (10 Rollers)	W.T.:0.8~1mm, I.D.:≤3.17mm	0.00011~32
BT100S-D3	2	DG6-2/DS6-2 (6 Rollers)	W.T.:0.8~1mm, I.D.:≤3.17mm	0.00016~49
BT100S-D4	2	DG10-2/DS10-2 (10 Rollers)	W.T.:0.8~1mm, I.D.:≤3.17mm	0.00011~32
BT100S-D5	4	DG6-4/DS6-4 (6 Rollers)	W.T.:0.8~1mm, I.D.:≤3.17mm	0.00016~49
BT100S-D6	4	DG10-4 /DS10-4 (10 Rollers)	W.T.:0.8~1mm, I.D.:≤3.17mm	0.00011~32
BT100S-D7	1	DT10-18	13#,14#, W.T.:0.8~1mm, I.D.:≤3.17mm	0.0023~64

BT100S-D8	2	DT10-28	13#,14#, W.T.:0.8~1mm, I.D.:≤3.17mm	0.0023~64
BT100S-D9	4	DT10-48	13#,14#, W.T.:0.8~1mm, I.D.:≤3.17mm	0.0023~64
BT100S-D10	1	YZ15/YZ15T	13#,14#,19#,16#,25#,17#,18#	0.005~641
BT100S-D11	1	YT25	114#,116#,15#,24#,35#,36#	0.024~750
BT100S-D12	2	2 x YZ15/YZ15T	13#,14#,19#,16#,25#,17#,18#	0.005~641
BT100S-D13	2	YT25 x 2	114#,116#,15#,24#,35#,36#	0.024~750
BT100S-D14	2	DT15-24	19#,16#,25#,17#	0.067~560
BT100S-D15	4	DT15-44	19#,16#,25#	0.067~330